KERN & SOHN GmbH
Ziegelei 1, 72336 Balingen-Frommern
Germany
Tel: +49 [0]7433 9933-199
service@kern-sohn.com

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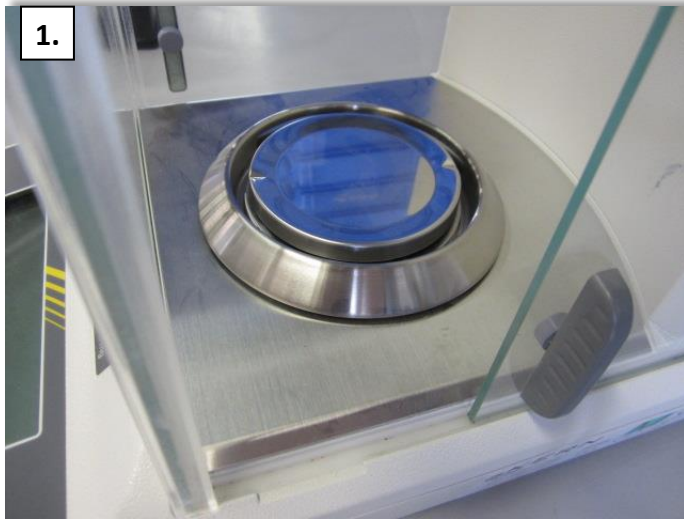
Important packaging instructions

Dear customer,

to prevent your measuring devices from damage during the transport, please consider the following packing instructions.

As measuring devices are always very delicate, please make sure that you use a covering box with enough cushioning material. It would be the best, if there is at least 15 cm - 20 cm cushioning material between the measuring device and the covering box to each side.

The following instructions apply especially for devices like analytical balances, moisture analyzers and precision balances.

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Remove **all** loose parts, such as weighing pan, pan support, protective ring, possible weighing space covers etc. from the weighing space.

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Pack the loose parts including the power supply unit in packing bags or bubble wrap and put them in a fitting box.

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Secure each sliding doors (top, on the right and left) against inadvertent opening with an adhesive tape

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Select a stable covering carton of a size that provides a crumple zone of at least 10 cm - 15 cm on all sides of the balance, that is, underneath, to the left and right, in front and behind and above the balance.

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Fill the covering carton base with a 10 cm - 15 cm layer of polystyrene chips or suitable spacing material (do not use air cushions or

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Place the balance **in a standing position** in the covering box on a bed of polystyrene chips or suitable spacing material (approx. 10 cm - 15 cm distance from the base).

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Place the small carton containing the loose parts (weighing pan, pan support, protective ring, power supply unit etc.) in between the back of the device and the wall of the covering box.

Use plenty of padding material all around the device, such as special packing paper or crumpled paper.

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Fill the covering box with padding material right up to the upper edge of the balance.

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Fill the remaining space right up to the covering box upper edge (approx. 10 cm - 15 cm) with polystyrene chips or some other suitable spacing material (do not use air cushions or solid, non-springy materials). Please enclose a delivery note including contact details and a detailed description of faults.

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